

HEMATOXYLIN EOSIN

1. Deparaffinization:
2. 6 min xylene 1
3. 6 min xylene 2
4. 6 min 100% ETOH
5. 6 min Xylene 2
6. 6 min 100% ETOH
7. 6 min 95% ETOH
8. 6 min 80% ETOH
9. 6 min 50% ETOH
10. Wash in distilled water
11. Incubate for 5 min in 1X TRIS or in PBS
12. Stain the nuclei with hematoxylin (DakoS2020) (3 min)
13. Wash in H₂O for at least 10 minutes to turn the violet color into blue
14. Quick wash in distilled H₂O
15. Color the cytoplasm with eosin (3 min)
16. Wash in distilled H₂O
17. Ethanol (96°) for 5 min (3 steps)
18. Xylol (Xylene)
19. Mounting with Canada balsam

Autor: M. Todaro

01/02/2015

Approved: Giorgio Stassi

